Center for Lasers and Photonics IIT Kanpur

Femtosecond laser laboratory

Precautions

- 1. It is mandatory to wear the lab footwear while entering the lab. Cell-phones should not be used inside the lab.
- 2. Equipment to be kept ON always
 - a. DT-506 main power switch
 - b. Power to ORC power supply
 - c. Water circulator chiller (THERMOTEK).
- 3. Room temperature should be maintained at $(18\pm2)^{0}$ C and humidity below 55%.
- 4. When the experiment is going on (or when the laser is kept ON), no more than 3 people can be present in the laser lab.
- 5. Goggles should be worn while working on the system.
- 6. Do not plug in or plug out any connections from the laser or any other device without prior permission.
- 7. Do not disturb any optical setup from optical table and/or with Micromachining system.
- 8. The assigned TA will change the output power of CPA; any other change in settings of DT 506 unit is strictly prohibited.
- 9. Be cautious always inside the lab. Keep the lab clean. Avoid any lengthy discussion with others while working inside the lab.
- 10. Ensure that all systems are shut down or stopped as per the procedure before leaving the lab.